

# MEASUREMENT OF DAIRY AND FOOD EMULSIONS USING LASER DIFFRACTION

## INTRODUCTION

The particle size of the fat droplets present in dairy and other food emulsions is important in defining properties such as flavour release, mouth feel and the emulsion stability. Large emulsion droplets can lead to poor flavour release, a greasy mouth feel and poor stability due to creaming. Emulsification to a smaller droplet size tends to reduce creaming and improve the taste of a product. However, in doing this a balance is required, as decreasing the particle size increases the available surface area, which in turn can lead to flocculation if the emulsifier concentration is not controlled.

In other products, such as ice cream, the particle size of the fat droplets is important in defining structural characteristics. Aggregated fat clusters are known to be involved in the stabilisation of the air cells within whippable dairy products. The formation of these clusters can only be achieved by the controlled destabilisation of the fat emulsion. Thus, a knowledge of the particle size is important in defining the functionality and taste of different food emulsion products.

## EMULSION MEASUREMENTS

The Malvern Mastersizer 2000 provides an excellent tool to the food scientist for the characterisation of food emulsions. Its wide dynamic range (0.02 2000 microns) permits both fine emulsion droplets and larger flocculated or coalesced droplets to be characterised. This range also allows for the measurement of large protein micelles, such as casein, enabling the interaction between the protein and emulsified fat phase to be understood.

## CHARACTERISING DIFFERENT MILK PRODUCTS

The particle size of milk products can be easily assessed using laser diffraction, allowing changes in the fat-phase to be detected. An example of this is shown in figure 1 where typical results for full-fat (3.6% fat), semi-skimmed (half-and-half, 1.7% fat) and skimmed milk (0.1% fat) are shown. As can be seen, two modes can be detected in each sample, one relating to the fat phase and one relating to free casein micelles. In moving from full fat to skimmed milk the relative proportions in each mode changes, tracking the reduction in fat content.

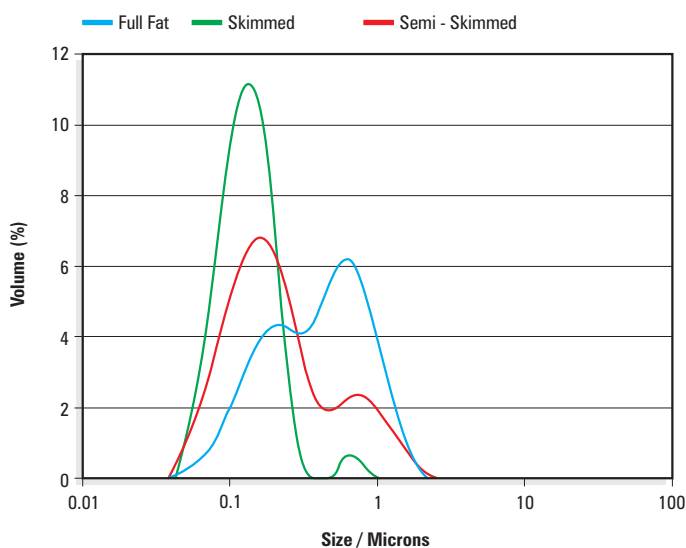


Figure 1: Size distributions recorded for Full Fat, Semi-Skimmed (Half-and-Half) and Skimmed Milk.

### MILK HOMOGENISATION

During processing, milk emulsions are normally homogenised in order to reduce creaming during storage. Laser diffraction can be used to track the progress of homogenisation, as shown in figure 2.

During the homogenisation of a milk emulsion (red curve, figure 2), a decrease in particle size is initially observed as the homogenisation pressure is increased. However, at high pressures the observed decrease becomes less pronounced. This is due to fat-cluster formation caused by bridging of the casein protein between the fat droplets within the emulsion. This occurs when the surface area of the fat droplets becomes too large to be covered by the available protein. Formation of these fat clusters can be inhibited using an appropriate “casein-dissolving” solution<sup>[1]</sup>. Adding this to the milk emulsion disperses the fat clusters, yielding a smaller particle size (blue curve, figure 2).

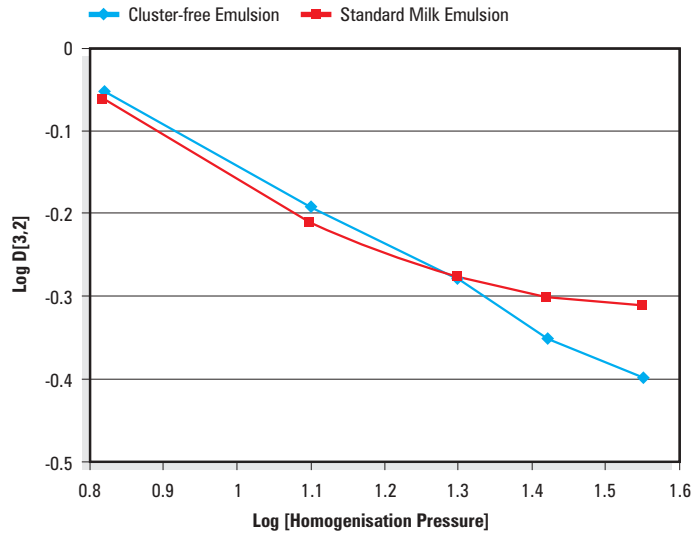


Figure 2: Variation of the D[3,2] with homogenisation pressure for a standard milk emulsion and a cluster-free emulsion containing the “casein-dissolving” solution<sup>[1]</sup>.

### EMULSION STORAGE

The behaviour of dairy emulsions during storage can also be related to particle size. Often emulsions such as cream liqueurs are found to increase in viscosity and even gel during prolonged storage<sup>[2]</sup>. Figure 3 shows how the Dv90 (particle size below which 90% of the volume of droplets exists) varies as a function of the viscosity measured over time for different liqueurs. Changes in the Dv90 can be used to detect the appearance of large particles. As can be seen, a direct correlation is observed between the Dv90 and the viscosity, with a move to the coarser particle sizes as the viscosity increases. This is caused by the formation of a flocculated droplet network.

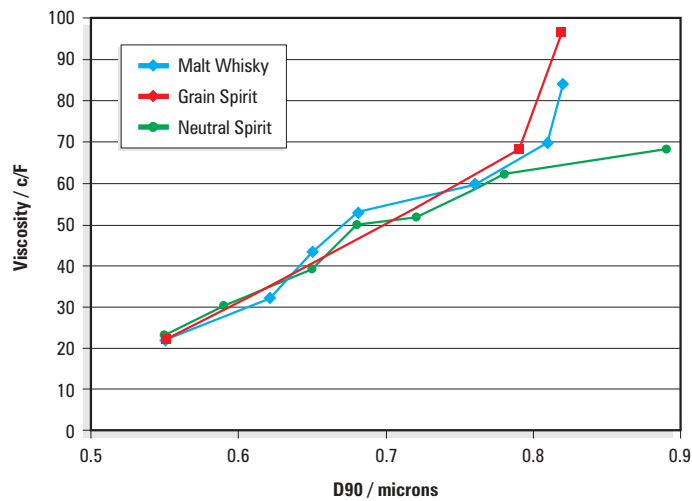


Figure 3: Variation in the particle size observed during the storage of cream liqueurs<sup>[2]</sup>.

## MILK POWDER RE-HYDRATION

Milk products are often spray dried before being shipped and reconstituted. The process of reconstitution of the spray-dried powder is an important factor in the production of many foodstuffs. This can be followed using laser diffraction.

Figure 4 shows the evolution of the particle size of an aqueous solution containing 5%w/v milk powder. The initial size of the powder was relatively large ( $> 10$  microns). Over time, samples of the solution were taken and measured. As can be seen, a mode at very fine particle sizes was observed as the larger powder mode decreased in volume. This fine mode relates to protein micelle formation during the re-hydration of the powder. Hydration is initially rapid but then slows down dramatically, with the process taking several hours to complete. In this case skimmed-milk powder was used, thus no fat was detected.

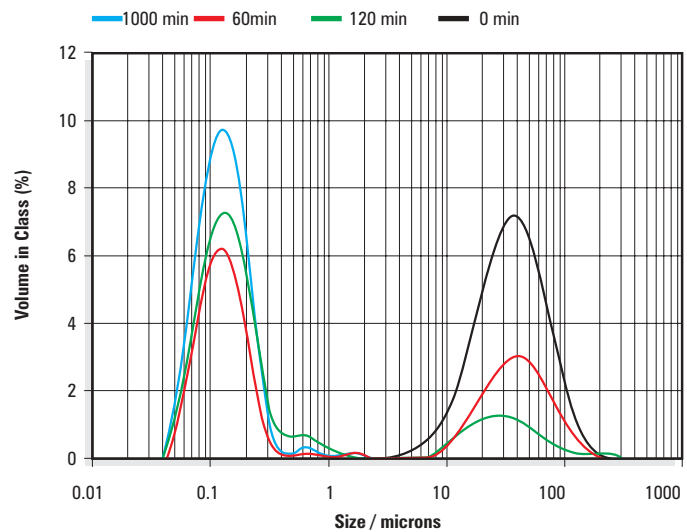


Figure 4: Milk powder reconstitution following using the Mastersizer 2000.

## CONCLUSIONS

The particle size of dairy and other food emulsions is an important factor in defining both structural and sensory characteristics. Measurements using the Mastersizer 2000 can be used to understand the changes in size that occur during production and storage of dairy products. This in turn can lead to a better understanding of how product formulation and performance are linked.

## ACKNOWLEDGEMENTS

Many thanks go to Catharina McCrae of the Hannah Research Institute for providing some of the data presented here.

## REFERENCES

- 1, McCrae, C.H. & LePoetre, A., "Characterisation of dairy emulsions by forward lobe laser light scattering - application to milk and cream", In. *Dairy Journal* (1996) 6: 247-256.
- 2, Muir, D.D., McCrae-Homsma, C.H. & Sweetsur, A.W.M1 "Characterisation of dairy emulsions by forward lobe laser light scattering - application to cream liqueurs", *Milchwissenschaft* (1991) 46: 691-694

**Malvern Instruments Limited**  
Enigma Business Park  
Groveswood Road  
MALVERN, Worcs.  
WR14 1XZ  
United Kingdom  
Tel: + 44 (0) 1684 892 456  
Fax: + 44 (0) 1684 892 789

**Malvern Instruments Inc.**  
10 Southville Road  
SOUTHBOROUGH  
MA 01772  
U S A  
Tel: + 1 (508) 480 0200  
Fax: + 1 (508) 460 0692

**Malvern Instruments Japan,**  
A Division of Spectris KK.  
South Bldg. 459,  
Kobe International  
Business Center (KIBC)  
5-5-2 Minatojima-minami-machi,  
Chuo-ku, Kobe-shi,  
Hyogo Prefecture, 650-0047, Japan  
Tel: + 81 (0) 078-306-3806  
Fax: + 81 (0) 078-306-3807

**Malvern Instruments GmbH**  
Rigipsstraße 19  
71083 HERRENBERG  
Germany  
Tel: + 49 (0) 70 32 97 77 0  
Fax: + 49 (0) 70 32 77 85 4

**Malvern Instruments SA**  
Parc Club de L'Université

30, Rue Jean Rostand  
91893 ORSAY CEDEX  
France  
Tel: + 33 (1) 69 35 18 00  
Fax: + 33 (1) 60 19 13 26

**Malvern Instruments Nordic AB**  
Box 15045  
Vallongatan 1  
750 15 UPPSALA  
Sweden  
Tel: + 46 (0) 18 55 24 55  
Fax: + 46 (0) 18 55 11 14

**WWW.MALVERN.CO.UK**